



Hutner's salts (from medium 590)

Nitrilotriacetic acid (NTA)	10.000	g
MgSO ₄ x 7 H ₂ O	29.700	g
CaCl ₂ x 2 H ₂ O	3.335	g
(NH ₄) ₆ Mo ₇ O ₂₄ x 4 H ₂ O	9.250	mg
FeSO ₄ x 7 H ₂ O	99.000	mg
"Metals 44"	50.000	ml
Distilled water	950.000	ml

1. Dissolve the nitrilotriacetic acid.
2. Adjust pH to 7.0 with KOH (~7.3 g).
3. Dissolve other salts separately and combine.
4. Adjust pH to 6.8 with NaOH or H₂SO₄.